manual

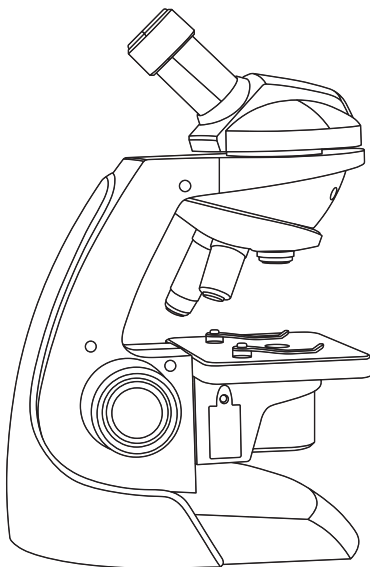
Microscope science kit 800x

Item: 47151

EN

NO

SV


www.playknowlogy.se
Box 50435 Malmö
Sweden
2022-09-27

Playknowlogy®

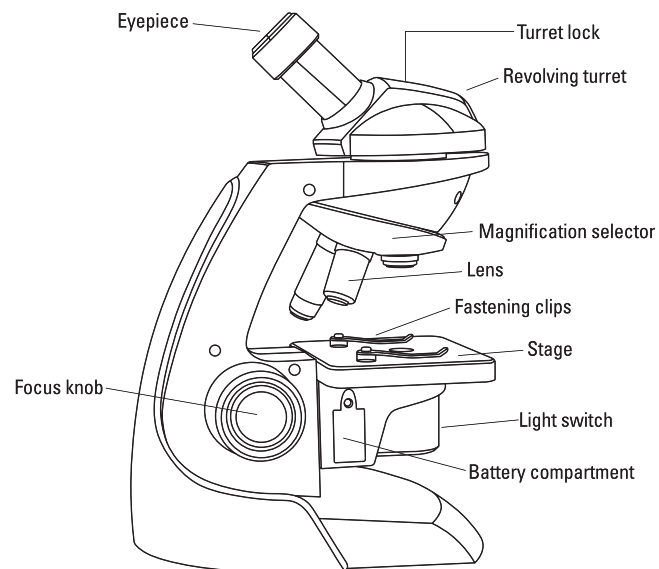Overview

Specifications

Magnification:	80x/200x/800x
Batteries:	3x LR44 (1.5 V)
Dimensions:	244x173x116 mm
Weight:	412 g

Included

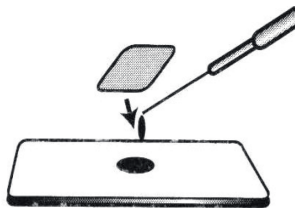
Microscope 80x/200x/800x, 12x blank labels, 12x cover glasses, 10x blank slides, 3x LR44 batteries, 2x collecting vials, 2x prepared slides, 1x pair of tweezers, 1x smartphone adapter, manual.

Use

The microscope is best used for samples that are thin and/or crystalline. If the sample is transparent, you can add drops of methylene blue, eosin or other dyeing solutions to the slide before placing the cover glass to better observe the sample. Samples of wool, pollen, salt or similar does not need a cover glass.

Preparing a temporary slide

1. Wipe the slide and cover glass clean with a soft paper towel. If needed, clean with a damp cloth to remove leftover dye or stubborn dirt.
2. Thin and flatten the sample using a scalpel or razor blade.
3. Place the sample as centered as possible on the blank slide using the tweezers.
4. Add a drop of water to the sample using a dissection needle or similar instrument. If the sample is transparent add your dye at this stage.
5. Cover the sample with the cover glass, make sure any pockets of air or bubbles are removed and remove any excess water or dyeing solutions with a paper towel.

Prepare a permanent slide

To prepare a permanent slide you can add a few drops of transparent adhesive glue or gum media around the sample during stage 4 in the above preparations. Leave the slide to dry for up to 24 hours before you handle the prepared slide, depending on the glue.

Using the microscope

1. Place the prepared slide on the stage and fasten it using the fastening clips.

2. Select the desired magnification strength by turning the magnification selector. You hear a click-sound when one of the 3 magnification settings is properly selected. Inspection of the sample is usually made at a lower setting, make sure the sample is centered below the lens and start increasing magnification.

3. Slide the light switch to  to turn on the light located under the stage.

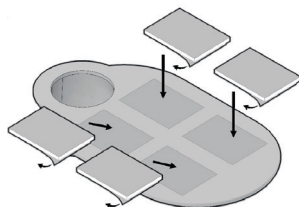
4. Turn the focus knob clockwise to move the lens as close as possible to the prepared slide. Look through the eyepiece and slowly turn the focusing knob counter-clockwise until the sample is in clear view and in focus. You can adjust the direction of the eyepiece by unlocking the turret lock and turning the revolving turret.


Attaching a smartphone

Instead of looking directly into the eyepiece you can use a smartphone with the included smartphone adapter.

1. Remove the protective film on one side of the four included adhesive tape pads. Attach them to the adapter.

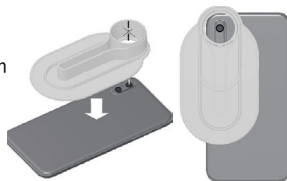2. Remove the top of the adhesive tape and align the center of the viewing tube to the smartphone camera. If misaligned you can lift the adapter gently and apply it again. Once aligned apply slight pressure to the adapter to secure it to the smartphone.

Note! The attachment area needs to be clean from fingerprints and dust, clean with a damp paper towel if needed and let the surface dry. Do not touch the adhesive surface. Where available a phone case can be attached with the adapter for further use.

3. Attach the viewing tube to the microscope eyepiece. Make sure the adapter is fully attached to avoid light leaking in for best picture result. Use the camera app on your device to view the sample through your microscope.

Replace batteries

If the stage lighting is weak, the batteries need to be replaced or inserted. Open the battery compartment with a screwdriver and insert 3x LR44 (1.5V) batteries. Mind the polarity. Close the battery compartment and fasten it with the screwdriver.

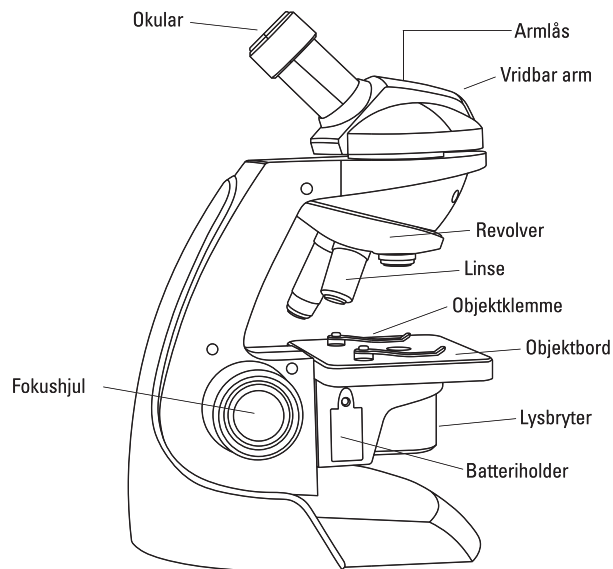
Safety information

Only adults should install and replace batteries.

Not suitable for use by children under 3 years as the product contains small parts that poses a choking hazard. Only for use under direct supervision of an adult. This toy contains sharp parts with sharp edges.

Remove the batteries if the microscope is not used for a long time.

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